

Settings > User

Account Preferences [Edit account preferences](#)

Username: [redacted]

Name: [redacted]

Email: [redacted]

Password: (hidden)

Institution:

Time zone: Pacific Time (US & Canada)

Research summary: [Enter one now](#) so colleagues in your CDD Vaults can learn more about activities.

Search across all vaults: false

Set time zone


Search across one or all vaults

Email Alerts

Type	Project	Frequency	Format	Protocols
New molecule	Any project	weekly	summary	
New batch	Any project	weekly	summary	
New protocol data	Any project	weekly	summary	Any protocol Preview

Alerts send email notifications when certain events occur.

Notification of additions to vault.
Note: No proprietary data is sent.

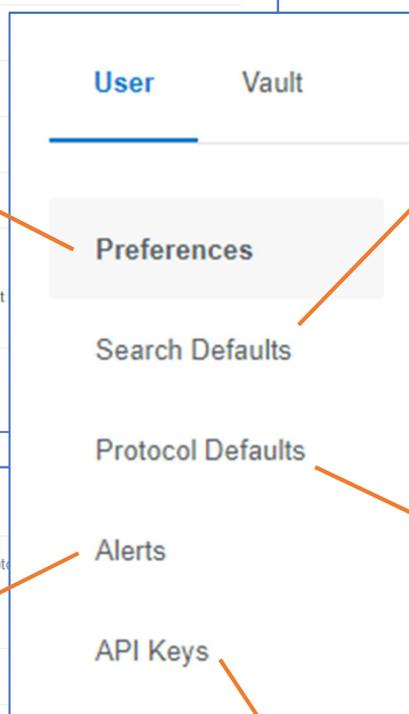

Search Defaults [Edit search defaults](#)

Detail level: Details

Displayed readouts: Matching

Dose-response plots scale: Per Run

Image size: Large

Column widths: 25 characters

Structure Fields: Structure, CDD Number

Batch Fields: Name, Salt, Scientist, Initial Amount, Purity, Comments, NMR

Specify options and batch fields to be displayed for every search

Protocol Defaults

[Add Protocol:](#) (select protocol)

Inhibition: All None

Run Date Inhibition (%) Avg Inhib (%)

Project Name SEM (%) Avg SEM

[Update protocol defaults](#) or [cancel](#)

Specify protocol fields to be displayed by default for each protocol

Your API Keys

Name	Role	Token
DemoRead	read only	[hidden]

[input field] read only [dropdown] [will be generated]

[Save](#) or [cancel](#)

After creation, token cannot be retrieved

Your new token named Delete is below. Please copy it to a safe place - you won't be able to retrieve it later.

[redacted token]

Note : you must protect this token as if it were a password.

Setting > Vault

Vault Settings

Vault name: Readout_definition

Primary contact: Readout: experiment

Principal investigator: Run: run

Institution: Molecule: compound

Website URL: Batch: lot

Next Registration Sequence Identifier: HA-0000515

Registration System Settings: Multiple compounds with the same structure are NOT allowed. To enable this feature (e.g. to archive structures with different but unique identifiers) please contact support@collaboratedrug.com.

Email alerts: summary only

Project association via data import: acceptance

External link setting: Ban external links Allow external links Allow external links and do not show a warning page

Session timeout: 120 minutes

Disable inactive members: Never

Authorized IP ranges:

Block CDD staff access: false

User Vault

Projects

Members

Terminology

General

Batch Fields

ELN Fields

3 Projects [Create a new project](#)

▶ CRO 1 [View data](#) [Delete](#)

▶ Project ABC [View data](#)

▶ Project XYZ

5 Members [Add a member](#)

Name	Access	Date Added	Projects	Email	API Keys	Actions
Heather Armaiz	vault admin	11/30/2017	3 Projects	heather@collaboratedrug.com	export, full access	edit delete
Maria Edwards	biology full access	6/5/2018	1 Project	harmaiz@collaboratedrug.com		edit delete
	read-add					edit delete
	full access	5/1/2018	2 Projects	whitney@collaboratedrug.com		edit delete

Batch Fields [Add/Edit Batch Fields](#)

Name	Data Type	Must be Unique	This Field
Batch Name			is required when creating batches
Date	Date		is optional
Scientist	Text		or Vendor is required
Notebook	Text		or Notebook is required
Vendor	Text		or Notebook is required
Sample ID	Text	✓	is optional
Initial Amount	Number		

Note: Uniqueness is compared across all batch fields within the vault.

ELN Fields [Add/Edit ELN Fields](#)

Name	Data Type	Must be Unique	Is Required
Project	Project		✓
External Journal ID	Text		
Experiment	Pick List		
MSDS	File		

Use your terminology in your Vault

Create projects and assign members to them

Read-Add or higher privileges needed

Create new members

Manage Vault members

Vault Admin privileges needed for all pages except Projects

General Vault setting

Create batch fields

Metadata fields for ELN entries

Specify batch field is:

- Unique value across Vault
- Required or optional