

Recommendations and tips

Modify terms used in CDD Vault

Settings > Vault > Terminology

Terminology [Edit terminology](#)

Vault Admin role needed

Suggested options for Protocol

- Results
- Data

Suggested options for Molecule

- Entity
- Construct
- Antibody
- Exosome
- Cell lines
- etc.

Entry:	entry
Protocol:	results
Readout_definition:	readout definition
Readout:	readout
Run:	run
Molecule:	entity
Batch:	batch

Help · Log out

Heather Arnaiz ▾

Entities Results Plates

Manually attach files

[Upload a file](#) · [Select an existing file](#)

Browse to a file

Associate a file that has been previously uploaded

Remove structure from search result table

Structure Fields · All · None

Structure

Uncheck Structure Fields

Setting > User > Search Defaults > Edit search defaults
OR
Click "Customize your report" from search result table

Best practices

Storing data in CDD Vault

Type of data	Entity	Entity descriptors	Result data	Procedures
	Antibody Construct Exosome Protein Cell line etc	Sequence Name PDB ID Isolation method Mutation etc	IC50 Western blot Validation Microscopy etc	Experiment details
Storage location in CDD Vault	<ul style="list-style-type: none"> Molecule <ul style="list-style-type: none"> Register without a structure Use Synonym to enforce uniqueness Recommend unique sequence, name or other identifier Multiple synonyms allowed Can be used to upload results 	<ul style="list-style-type: none"> Batch fields <ul style="list-style-type: none"> Create batch fields to store the information <p style="text-align: center;">OR</p> <ul style="list-style-type: none"> Protocol <ul style="list-style-type: none"> Create one or more protocols to store the information <p>Protocol = container to store data and perform calculations</p>	<ul style="list-style-type: none"> Protocol <ul style="list-style-type: none"> Create one or more protocols to store the results and perform desired calculations. <p>Protocols are very flexible. A protocol can be created for one readout or many.</p>	<ul style="list-style-type: none"> ELN <ul style="list-style-type: none"> Store details of experiment Link to related ELN entries Link to results and entities